

PRODUCT DATA SHEET

Bioworld Technology CO., Ltd.



DAB Chromogen Kit

Cat No.: BD5038

Introduction

This product is prepared for Immunohistochemistry, Immunoblot or Dot Blot. The positive cells will be an insoluble brown sediment. The insoluble brown sediment in form of bands or dots.

Reagents

A: PBS	60ml
B: Substrate Reagent	3ml
C: DAB Chromogen	3ml

Application

Used for IHC and WB Color reaction.

Storage & Shelf life

Store at 2-8°C for short time; -20°C for long time. Each component is stable for up to 12 Months.

Procedure

1. In order add: 1ml PBS (Vial A), 50ul Substrate Reagent (Vial B), 50ul DAB Chromogen (Vial C), mixing vial shortly before incubation;
2. Apply 50-100ul DAB Substrate Reagent to each slide, incubate about 3-5 minutes;
3. When sufficient staining has been achieved rinse slides in PBS for 3 minutes for 3 times.

RESEARCH USE

For research use only, not for use in diagnostic procedures.

MADE IN CHINA

Bioworld Technology, Inc.

1660 South Highway 100, Suite 500 St. Louis Park, MN
55416, USA.

Email: info@bioworld.com

Tel: 6123263284

Fax: 6122933841

Bioworld technology, co, Ltd.

No 9, weidi road Qixia District Nanjing, 210046,
P, R.China.

Email: info@biogot.com

Tel: 0086-025-68037686 Fax: 0086-025-68035151